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
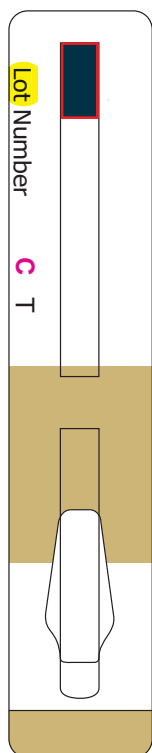
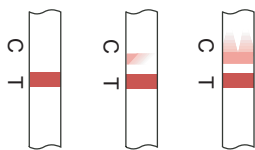
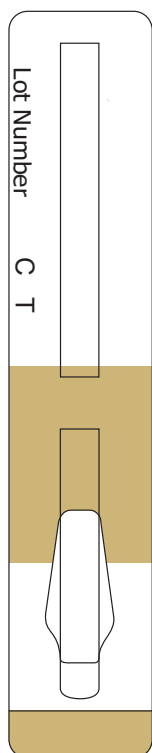
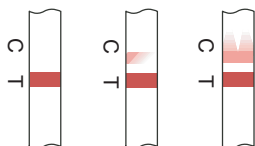
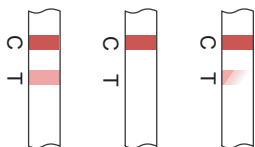
POSITIVE


NEGATIVE

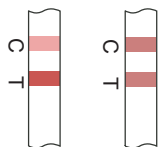Note: Visual interpretation is
NOT recommended.

PROCEDURE

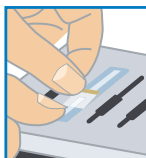①

- Verify incubator temperature is $56 \pm 1^{\circ}\text{C}$.
- Shake milk sample and label test strip.
- Place strip in incubator and use tab to peel tape back to "Peel to here" line.

②

- Slowly pipet 300 μ l sample into sample compartment at ROSA® Incubator indicator line (as shown).
- Reseal tape over sample compartment.
- Close lid on incubator and tighten latch.
- Incubate for 3 minutes, but not more than 4 minutes.

③

- Remove tests from incubator.
- Visually inspect C (control) line for even development.
- Read with ROSA® Reader.
- Insert clean and visually valid test strip in ROSA Reader and read results on MRL channel in 2-line mode (slow-blinking).

© 2012 Charm Sciences, Inc. Charm, ROSA names are registered trademarks and test strip product packaging is trade dress, of Charm Sciences, Inc. MRK

SL® Level

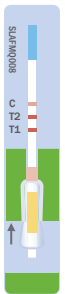
Charm® 3 SL3
3-Minute Beta-lactam


SL
Beta-lactam


SLAFMQ
SL Aflatoxin Quantitative

FLUSLBL
Beta-lactam / Flunixin

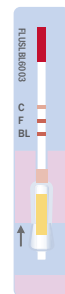
SLAFM
SL Aflatoxin Visual Test

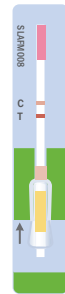
SULF
Sulfonamides

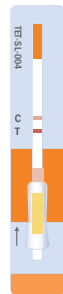
TET-SL
Tetracycline

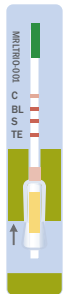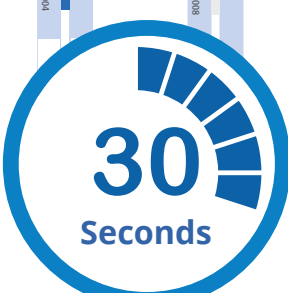TRIO
Beta-lactams,
Sulfonamides, and
Tetracyclines


**30
Seconds**

SL / MRL Levels

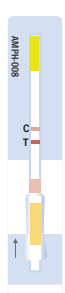
AMPH
Amphenicol


TET
Tetracycline

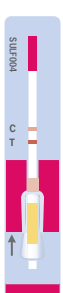
SULF-MRL
Sulfonamides


QUIN
Quinolones

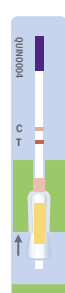
SL-KIWI
Beta-lactam

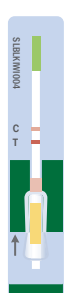
CAP
Chloramphenicol

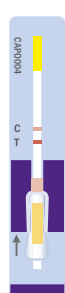
STREP
Streptomycin

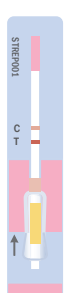NEO-STREP
Neomycin/
Streptomycin

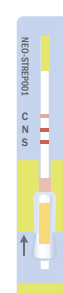PIRL
Pirlimycin

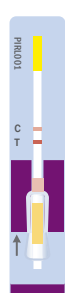
MRL Level


AUSBL
Beta-lactam

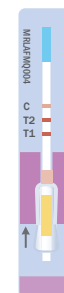
MRLAFMQ
MRL Aflatoxin
Quantitative

MRLBL3
Beta-lactam
3-Minute

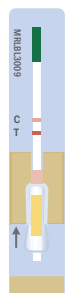MRLBL
Beta-lactam

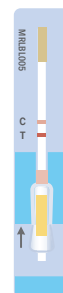MRLBL1
Beta-lactam
1-Minute

MRLBLTET
Beta-lactam
Tetracycline

MRLBLTET2
Beta-lactam
Tetracycline
2-Minute


MRLBLRFTET2
Beta-lactam
Tetracycline
2-Minute


PLABLTET
Beta-lactam
Tetracycline
8-Minute

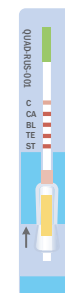QUAD
Chloramphenicol,
Beta-lactams,
Tetracyclines,
and Streptomycin

QUAD 1
Sulfonamides
Beta-lactams,
Tetracycline,
and Quinolones

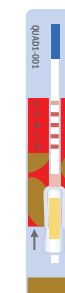QUAD2
Lincomycin/Pirlimycin,
Tylosin/Tilmicosin,
Spiramycin, and
Erythromycin

QUAD3
Spectinomycin,
Kanamycin,
Neomycin, and Strep/
Dihydrostreptomycin

Charm Sciences, Inc.

659 Andover Street, Lawrence, MA 01843-1032, USA

Tel: +1.978.687.9200 | Fax: +1.978.687.9216 | Email: info@charm.com | www.charm.com

© 2017 Charm Sciences, Inc. Charm, SL, ROSA and Running Cow logo are registered trademarks, and test strip product packaging is trade dress, of Charm Sciences, Inc.
 MRK-170-008 Apr-2017